# RZR 1

**Betriebsanleitung**  
**Instruction Manual**  
**Mode d'Emploi**  
**Instrucciones de**  
**Empleo**

D	DEUTSCH	Seite	3 – 12
E	ENGLISH	page	13 – 22
F	FRANCAISE	page	23 – 32
ES	ESPAGNOL	pàgina	33 – 42

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#### **Niederspannungs-Richtlinie (73/23/EWG):**

EN 61010-1  
 EN 61010-2-051

Thank you for purchasing a Heidolph Instruments product. This item has been designed, made and inspected in compliance with DIN EN ISO 61010 for long-term and flawless operation.

## SUMMARY

<b>SUMMARY.....</b>	<b>13</b>
<b>STANDARD HARDWARE &amp; OPTIONS.....</b>	<b>14</b>
<b>GENERAL INFORMATION.....</b>	<b>15</b>
<b>SAFETY INFORMATION.....</b>	<b>15</b>
<b>ASSEMBLY.....</b>	<b>16</b>
<b>OPERATION AND CONTROLS.....</b>	<b>18</b>
<b>CLEANING &amp; SERVICING.....</b>	<b>19</b>
<b>DISASSEMBLE &amp; STORAGE.....</b>	<b>19</b>
<b>DISPOSAL.....</b>	<b>20</b>
<b>TROUBLESHOOTING.....</b>	<b>20</b>
<b>SPECIFICATIONS.....</b>	<b>20</b>
<b>WARRANTY, LIABILITY &amp; COPYRIGHT.....</b>	<b>21</b>
<b>QUESTIONS / REPAIR WORK.....</b>	<b>21</b>
<b>CE-DECLARATION OF CONFORMITY.....</b>	<b>22</b>

**Important information**

**Advice about power cord / mains supply**

**Caution: mandatory action**

**Caution: fire- and explosion hazard**

**Advice about maintenance / repair**

## STANDARD HARDWARE & OPTIONS

### Standard Hardware

	<b>Product</b>	<b>quantity</b>	<b>P/N 230V 50Hz</b>	<b>P/N 115V 60Hz</b>
	RZR 1	1	501-11000-00	501-11000-04
	support rod	1	22-02-14-01-14	22-02-14-01-14
	allen key	1	02-07-01-01-01	02-07-01-01-01
	chuck	1	509-01000-00	509-01000-00
	T-handle (for chuck)	1	02-07-03-01-01	02-07-03-01-01
	Instruction Manual	1	01-005-001-16	01-005-001-16

### Accessories & Stirring tools (optional)

	<b>Item</b>	<b>1st.1.1 P/N</b>
	stand	570-12000-00
	clamp	570-22000-00
	flexible stirrer coupling shaft	509-03000-00
	stirring paddle BR 10	509-10000-00
	stirring paddle BR 11	509-11000-00
	stirring paddle BR 12	509-12000-00
	stirring paddle BR 13	509-13000-00
	stirring paddle BR 14	509-14000-00
	prop-type stirring paddle PR 30	509-30000-00
	prop-type stirring paddle PR 31	509-31000-00
	prop-type stirring paddle PR 32	509-32000-00
	prop-type stirring paddle PR 33	509-33000-00
	turbine TR 20	509-20000-00
	turbine TR 21	509-21000-00
	half moon HR 18	509-18000-00
	anchor AR 19	509-19000-00
	prop-type stirring paddle PR 39	509-39000-00
	enclosure NS 29/32	509-09000-00

## GENERAL INFORMATION

-  Unpack your item carefully.  
Inspect for damage and report such damage or missing parts to your supplier right away.
-  Read your Instruction Manual carefully. Take time to save time when working with your product. Make sure that every user has read and understood the Instruction Manual.
-  Please store the Instruction Manual in a place easily accessible to every user.

IF ALL ELSE FAILS, READ THESE INSTRUCTIONS !

-  A so-called EURO-plug (DIN 49441 CEE 7/VII 10/ 16 A 250 V) is standard on all of the products.  
For the Continental US they feature a US-standard plug (NEMA Pub.No.WDI.1961 ASA C 73.1 . 1961 page 8 15A 125V).
-  For using the item in a country with deviating outlet / plug systems, we recommend to use approved adapters or to have an electrician replace the standard plug with one suiting your needs.
-  As shipped, the item features a protective ground wire. When replacing the original plug, make sure to reconnect this protective ground wire in the new plug !

## SAFETY INFORMATION

-  Please comply with all safety and accident-prevention regulations as in force for laboratory work!
-  Use extra care when working with flammable substances; refer to safety data sheets.
-  Use extra care when working in the vicinity of flammable and explosive substances. Motors are of non-sparking type, the item itself however is not explosion-proof.
-  When connecting your item with your local power supply, please make sure your item is designed for your local voltage; go by the data plate on the item.
-  Please connect your item with a protective-ground outlet only.
-  Turn your power switch OFF whenever the item is not used, or before disconnecting the plug.
-  Repair work is limited to technicians approved by Heidolph Instruments.
-  Your item needs a solid stand.

## ASSEMBLY

The RZR 1 comes with a support rod (B) a 4mm Allen key, chuck, chuck key and 2mm Allen key. The clamp, stand, and paddle have to be ordered separately. See accessories.

### Install stand

Loosen the Allen set screw (A) and insert the support rod (B) into the opening at the back of the stirrer (as in diagramm). Tighten set screw (a) with the Allen key provided.

### Install chuck

Place the chuck (D) over the hollow shaft of the stirrer and with a 2 mm Allen key tighten the Allen set screw (C) in the collar of the chuck. Ensure the chuck is safely fastened without damaging the hollow shaft.

Place desired paddle in chuck and tighten using the chuck key.


### Install stirrer on stand

Slide the support rod (B) with the stirrer into the opening of the clamp, adjust to desired position and tighten.


#### Attention:

- Ensure that the chuck is safely fastened
- Please strictly comply with current safety regulations for machines exposing rotating shafts.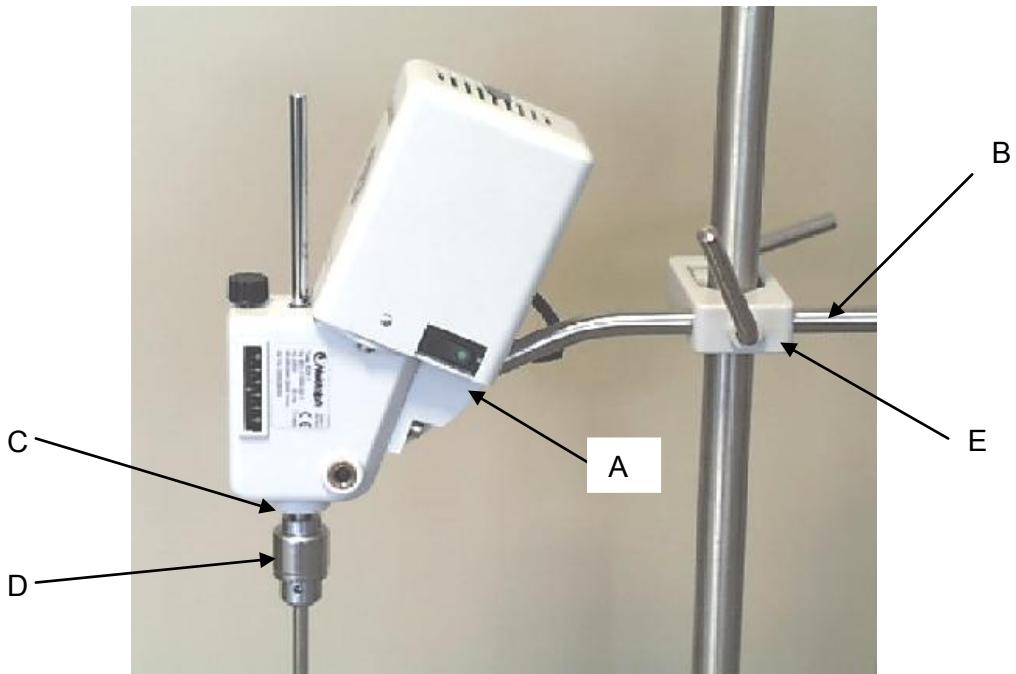


## Install stirring tool

Introduce stirring tool (N) into the chuck, put in place and secure with T-handle (M). Immersion depth of the stirring tools is adjusted by displacing the tool in and out in the open chuck (D).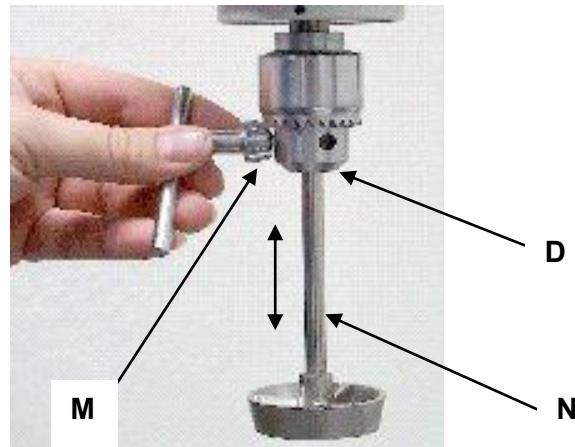

## Install flexible coupling shaft

Hold the flexible coupling shaft (O) in the chuck (D) by its metallic stud (P) and secure with T-handle (M). At this point introduce stirring tools (N) shaft end in collet (Q) and pull tight (cw)

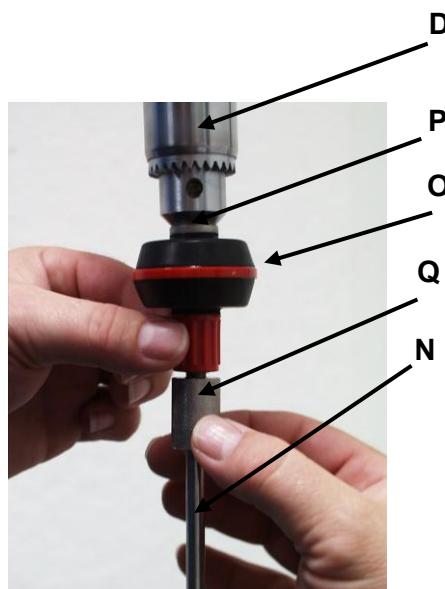## OPERATION AND CONTROLS

Before connecting power cord with mains outlet, make sure that:

- your item is designed for your local voltage and frequency (data plate on item).
- master switch is set to "0".

### Initial operation


Power supply is done by cable with plug with a length of 2 m. A rocker switch, located on the right side of the instrument, turns the stirrer off and on. A pilot light will indicate when the stirrer is on. The green pilot lamp within the ON/OFF switch should light up when the instrument is operational.

### Adjustment of speed


The RZR 1 has two speed ranges and two hollow shafts for mounting paddles depending on the desired speed and torque. In the upright position, the speed range is 280-2200 rpm. For the second speed range, the stirrer must be pivoted 90°. To do this, loosen the clamp holding, the support rod (e) and turn the stirrer. This shaft (1) has the speed range of 35-250 rpm. Adjustment within these two ranges is done with the black knob (F) at the front of the stirrer. For control purposes you find an indicating scale (G) at the side of the gear housing.

Shaft 1 : 35-250 min<sup>-1</sup>

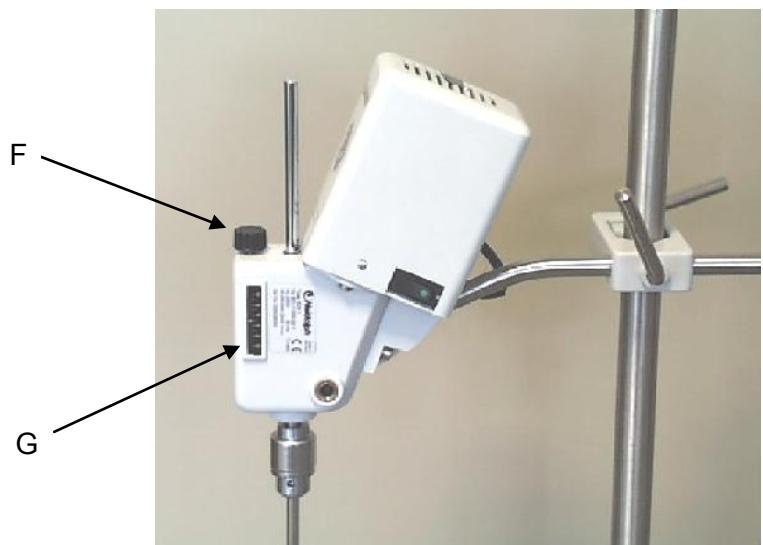Shaft 2 : 280-2200 min<sup>-1</sup>
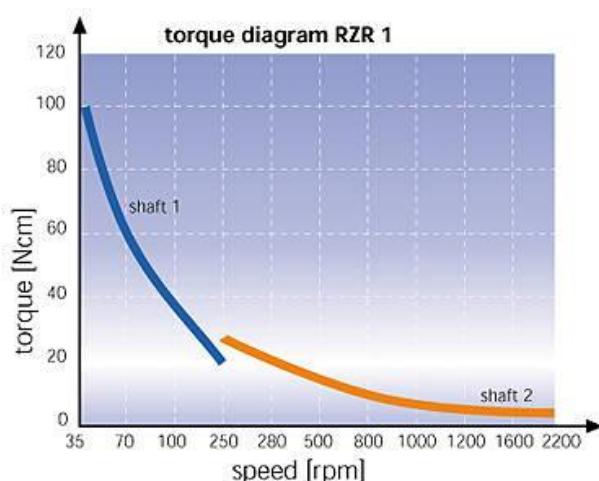
## CLEANING & SERVICING

**Cleaning:** wipe housing clean with a damp cloth (add some sort of mild liquid soap).

### Note

To avoid damage to the surface finish, avoid using chlorine bleach, chlorine-based detergents, abrasive substances, ammonia, rags or cleaning agents containing metal particles.

### Servicing

Motor and gear are fitted with attendance-free ball bearings and are constructed for continuous running. Subject to wear in such a gear is the friction lining, here the friction rubber. If the friction rubber has to be replaced one day, it can be done directly in the laboratory without considerable display. For this purpose the gear housing is opened by loosening of the fastening screws, and the main shaft with the friction wheel is removed. By means of a screw-driver the friction rubber can be taken away and replaced by a new one. Please notice that friction gears must not be lubricated or greased. When replacing the friction rubber you have therefore to pay attention that the friction cone must not be covered by an oil or grease film. For scouring you can use petroleum ether or spirit of wine. The construction of the gear is done in reversed sequence. Service life of the friction rubber depends on the precise pressure. The pressure of the friction cone acting against the friction wheel is adjusted correctly, if at adjustment-position "10" the motor comes to standstill at braking of the driving shaft. At letting loose of the driving shaft the motor must start immediately. This can be ascertained by taking a look through the slots of the motor hood.

The item is maintenance-free. Repair work is limited to technicians approved or appointed by HEIDOLPH Instruments. Please call your local HEIDOLPH Instruments Dealer or a HEIDOLPH Instruments Field Representative.

## DISASSEMBLE & STORAGE

### Disassemble

1. Turn the item OFF and disconnect mains plug.
2. Remove all of the hardware arranged around the item to ease disassembly.
3. Open chuck with T-handle and remove stirring tool
4. Unlock clamp and remove stirrer head from the head clamp
5. With the Allen key loosen threaded pin and remove chuck.

### Storage

1. We recommend to store the item and its components in its original box, or a similar container that offers adequate protection against damage in transit. Tape the box securely.
2. Store the item in a dry place.

### Caution

Do not jolt or shake the item during transport.

## DISPOSAL

For disposal, please comply with your local or national regulations.

Split by metal, plastic, etc.

Packing material to be treated as described above (material split).

## TROUBLESHOOTING

Work on electric and electronic components is limited to qualified personnel.

**Master switch on stirrer won't light**

1. check power cord and power supply

**Item won't work (master switch lighted)**

2. Thermal motor circuit breaker triggered by motor overload

Remedy:

Wait about 20 minutes, decrease load applied on shaker plate.

**Motor running, output shaft static**

3. rubber ring or friction cone are worn out

Remedy:

replace rubber ring or friction cone and adjust press load

## SPECIFICATIONS

**RZR 1**

Model	RZR 1
Motor power- input/output (W)	77/18
Number of gear stages	2
Speed range (1/min)	35-250 and 280-2200
Speed control	mechanically stepless, scale with guide numbers
Max.torque (Ncm)	100
Torque at 1000 min <sup>-1</sup> (Ncm)	8
Dimensions of unit (BxHxT) (mm)	71 wide; 250 high; 172 deep
Dimensions of stay bar (ØxL) (mm)	13x300
Weight (kg)	2.5
Ø Dia. of stirrer shaft (plug in) (mm)	max.8,0
Operating temperature range (°C)	0-40
Max.permissible relative humidity	max. 80% (avoid reaching dewpoint)
protection class	IP 20
Interference suppression standard	VDE 0875
Motor protection feature	overheating protection, self-resetting
Permissible operating time period	100%

## WARRANTY, LIABILITY & COPYRIGHT

### **Warranty**

Heidolph Instruments warrants that the present product shall be free from defects in material (except wear parts) and workmanship for three years from the date shipped off the manufacturer's warehouse.

Transit damage is excluded from this warranty.

To file for such warranty service, contact Heidolph Instruments (phone: +49-9122-9920-68) or your local Heidolph Instruments Dealer. If defects in material or workmanship are found, your item will be repaired or replaced at no charge.

Misuse, abuse, neglect or improper installation are not covered by this warranty.

Alterations to the present warranty need Heidolph Instruments' consent in writing.

### **Exclusion Clause**

Heidolph Instruments cannot be held liable for damage from improper use or misuse. Remedy for consequential damage is excluded.

### **Copyright**

Copyright in pictures and wording of the present Instruction Manual is held by Heidolph Instruments.

## QUESTIONS / REPAIR WORK

If any **aspect** of installation, operation or maintenance remains unanswered in the present Manual, please contact the following address:

For repair services please call Heidolph Instruments (phone: +49 - 9122 - 9920-68) or your local, authorized Heidolph Instruments Dealer.

### **Note**

You will receive approval for sending your defective item to the following address:

**Heidolph Instruments GmbH & Co. KG**  
**Technical Service**  
**Walpersdorfer Str. 12**  
**D-91126 Schwabach / Germany**  
**Tel.: +49 – 9122 - 9920-74**  
**Fax: +49 – 9122 - 9920-84**  
**E-Mail: [Service@Heidolph.de](mailto:Service@Heidolph.de)**

### **Note**

If you are based in the United States of America, please contact Heidolph US:

**Heidolph Instruments**  
**North America**  
**Phone: 1-866-650-9604**  
**E-Mail: [service@heidolph.com](mailto:service@heidolph.com)**  
**[www.heidolphNA.com](http://www.heidolphNA.com)**

### Safety Information

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact substance
- take proper protective measure to ensure the safety of our receiving and service personnel
- mark the pack IAW Hazardous Materials Act

### CE-DECLARATION OF CONFORMITY

We herewith declare that the present product complies with the following standards and harmonized documents:

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